



INSTALLATION INSTRUCTIONS

1. Switching on your unit

Once your unit arrives by freight, please wait for approximately 1 hour before switching on your unit to allow the oils in the compressor to settle after shipment. It is best to use a dedicated power source clearly marked "do not switch off. Once your unit has time to settle, plug in and turn on your unit.

2. Alarms

When the refrigerator is first plugged in, it will initially display a flashing "PF" alert on the Min-Max temperature display located on the left hand side of your unit informing the operator there has been a loss of power since its original set up in the manufacturing warehouse. To acknowledge this alert, press the red set button on the left Min-Max temperature display as per the image below. The refrigerator may take time to reach the desired temperature range, resulting in an additional alarm condition ("HI"). To silence the "HI" alarm, press the red set/mute button on the right hand electronic controller.



3. Register your cloud monitoring

Your Medisafe Plus unit comes complete with 12 months access to the Medisafe Plus Cloud monitoring system. As soon as your refrigerator is plugged in to a power source it will begin to record data automatically. To access this data please register your monitoring system on the Medisafe Plus Cloud Monitoring system portal.


Please complete your activation form by visiting www.medisafeplus.com.au/registration.pdf

www.medisafeplus.com.au

Nationwide Phone: 1 800 658 120

5. Equipment Serial Number

When registering your cloud monitoring system, please have your Equipment serial number handy. You can find your serial number on the white Avem Quirks sticker inside your fridge

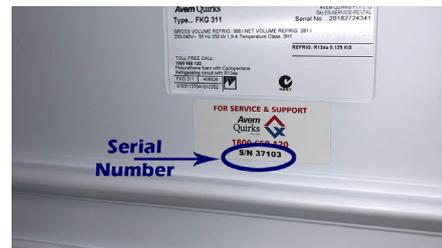

6. User Manual

A full user manual for your Medisafe Plus unit can be found by visiting

www.medisafeplus.com.au/usermanual.pdf

The user manual will explain how to read your data and download or print your temperature reports.

7. Service Support

It is recommended that your Medisafe Plus unit be serviced with an annual maintenance and calibration service each year. To book your service by an authorised technician please contact our office.

Each Medisafe Plus unit comes standard with manufacturers 24 months parts and labour warranty.

www.avemquirks.com.au/services/

E: Medisafeplus@avemquirks.com.au

P: 1 800 568 120

Supplied by

team medical
supplies

✉ P.O. Box 6245 Dural DC, NSW 2158

☎ 1300 22 44 50

📠 1300 22 44 60

@ info@teammed.com.au

🌐 www.teammed.com.au